



KOJAIR[®]

Biowizard Platinum SF

Biowizard Platinum MF

Biowizard Platinum DEF

Biowizard Platinum XL MF

Cleanwizard Platinum CW-V

Platinum products

Accessories

TABLE OF CONTENT

1	INSTALLATION OF EXHAUST DUCT CONNECTOR	3
2	ACCESSORIES LIST	5

1 INSTALLATION OF EXHAUST DUCT CONNECTOR


When working with cytotoxics or with other substances causing serious health hazards, the cabinet should always be connected to the exhaust duct. An exhaust duct connector is available as an optional accessory.

If the cabinet is connected to an exhaust duct, there has to be a separate exhaust fan in the duct. The cabinet does not protect the operator if the air, which is emitted, encounters resistance in the exhaust duct.

	NOTE! Use only the manufacturer's exhaust duct connector. The wrong type of exhaust duct connector might decrease the protection level of the cabinet. The manufacturer is not responsible for the function of the cabinet, if any other part is used.
--	---

The accessories needed for the exhaust connection are not included in the standard delivery. Accessories may be ordered with the cabinet as options or later on as accessories.

When cabinet is connected to external exhaust channel the exhaust air volume in delivery line must be 10 -15% higher than the cabinet's exhaust air volume. See section technical data.

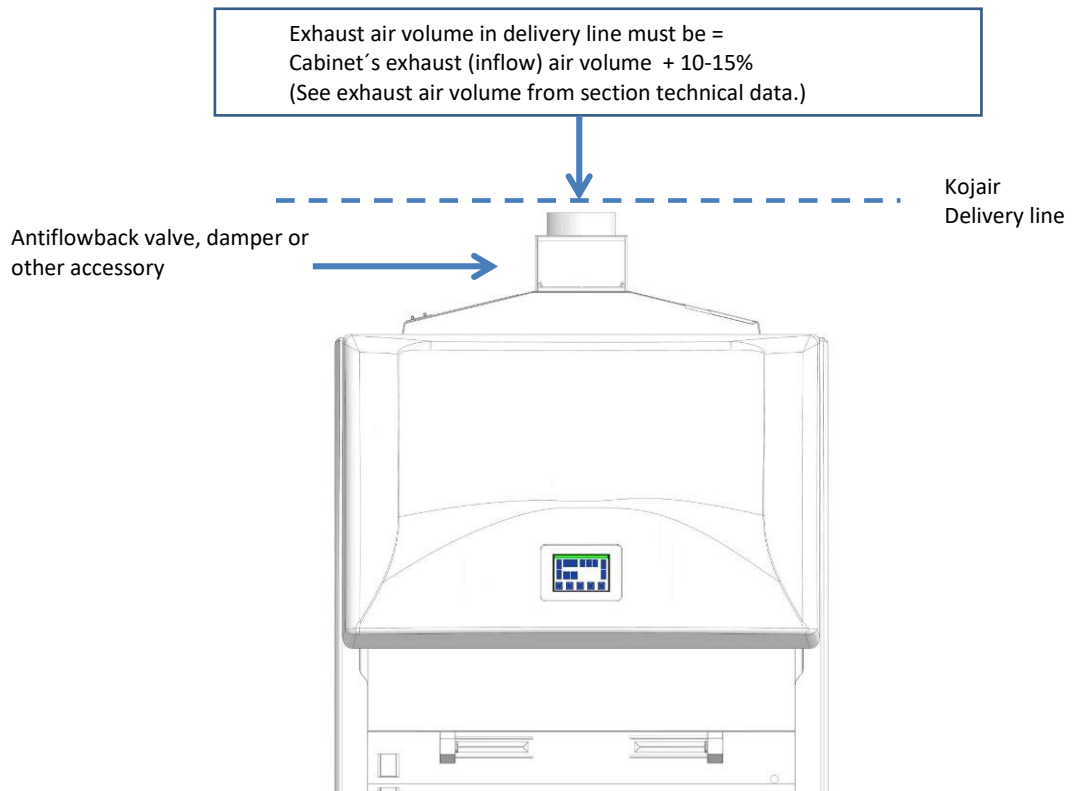


	NOTE! If new accessories are installed after installation tests, in some cases installation test have to carried out again. For more information please contact your local Kojair dealer.
--	--



NOTE! When connecting the cabinet to the exhaust duct an anti-flow back valve must be installed, because pressure changes in the exhaust duct might temporarily decrease the protection level. Installation is required by the EN 12469-norm.

If the anti-flow back valve is not ordered to your device, the anti-flow back valve must be located in the exhaust ventilation of the building.

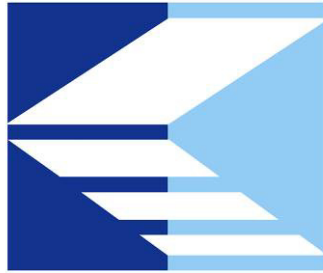

KOJAIR[®]

Kojair Tech Oy

Teollisuustie 3

35700 Vilppula

Finland

Switchboard: +358 3 471 7000

Fax +358 3 471 7077

info@kojair.com

www.kojair.com

